



**KERN & Sohn GmbH**

Ziegelei 1

D-72336 Balingen

E-Mail: [info@kern-sohn.com](mailto:info@kern-sohn.com)

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: [www.kern-sohn.com](http://www.kern-sohn.com)

# Installation Instructions

## Rechargeable battery

### **KERN FOB\_A07-A08**

Version 1.0

12/2010

GB

FOB\_A07-A08-IA-e-1010



# KERN FOB\_A07-A08

Version 1.0 12/2010

## Installation Instructions

### Rechargeable battery

Rechargeable battery	Platform
FOB-A07	180 x 170 mm
FOB-A08	215 x 215 mm (L)

## 1 Technical data

Rechargeable battery	Operating period	Loading time
FOB-A07	approx. 24 h	approx. 8 h
FOB-A08	approx. 30 h	approx. 14 h

## 2 Scope of delivery

Rechargeable battery	
FOB-A07	Battery block, 7.2 V; 1200 mAh
FOB-A08	Battery block, 7.2 V; 2000 mAh

## 3 To install rechargeable battery:

- ⇒ Disconnect the balance from the mains
- ⇒ Remove weighing plate

⇒ On the lower side of the balance loosen screws of the rechargeable battery compartment cover



⇒ Remove the rechargeable battery compartment cover



⇒ Plug connection of the battery block into connection of the balance





- ⇒ Insert the rechargeable battery block into the rechargeable battery compartment
- ⇒ Ensure that the cable is not squeezed.



- ⇒ Screw down the rechargeable battery compartment cover



- ⇒ Re-install the weighing plate

#### 4 Initial Commissioning

The rechargeable battery is loaded via the mains adapter of the balance. Before the first use, the rechargeable battery should be charged for 14 hours. Operating period of the rechargeable battery see table chap. 1 – Technical data.